# FUJIFILM

## Irvine Scientific

### Colcemid Solution

Catalog No. 9311

12 x 10 mL Kits

- 10 µg/mL in Hank's Balanced Salt Solution
- 10 µg/ml in Hank's Ausgeglichenere Salzlösung
- 10 µg/ml in Soluzione di Hank di Sale Equilibrato
- 10 µg/ml en Solución Salina Equilibrada de Hank's
- 10 µg/ml dans la Solution Sa line Équilibrée de Hank
- 10 µg/ml em Solução Salina Equilibrada Hank's
- 10 µg/mL σε διάλυμα αλάτων Hank (Hanks' Balanced Salt Solution/HBSS)
- 10 µg/ml v Hankově vyváženém solném roztoku
- 10 µg/ml i Hank's afbalancerede saltopløsning
- 10 µg/ml Hank's Balanced Salt Solution -suolaliuoksessa
- 10 µg/ml Henka līdzsvarotā sāls šķīdumā
- 10 µg/ml in Hanks gebalancēerde zoutoplossing
- 10 µg/ml w zbalansowanym roztworze soli Hanka
- 10 µg/ml în Soluție Salină Echilibrată Hank
- 10 µg/ml i Hanks balanserade saltlösning
- 10 µg/ml Hanki tasakaalustatud soolalahuses
- 10 µg/ml Hank-féle kiegyenlített sóoldatban
- Henko subalansuotas druskų tirpalas, 10 µg/ml
- Hanks Dengeli Tuz Solüsyonunda 10µg/ml
- 10 µg/ml v Hankovom vyváženom soľnom roztoku
- 10 µg/ml в балансиран солев разтвор на Ханк
- 10 µg/ml u Hankovoj izbalansiranoj otopini
- 10 µg/mL f'Soluzzjoni Salina Bilanċjata ta' Hank
- 10 µg/ml v Hankovi uravnoteženi solni raztopini

#### Glossary of Symbols\*:

	Catalog Number
	Lot Number
	Sterilized using aseptic processing techniques (filtration)
	Expiration: Year - Month - Day
	Caution, consult accompanying documents
	Consult instructions for use
	Storage Temperature 2-8°C
	Manufacturer
	CE Mark
	Emergo Europe - Prinsessegracht 20 2514 AP The Hague The Netherlands

\*Symbol Reference - EN ISO 15223-1, Medical devices – Symbols to be used with medical device labels, labeling.

#### REFERENCES

- L.S. Goodman and A. Gilman. *The Pharmacological Basis of Therapeutics*. 5<sup>th</sup> edition. p. 351. (Mac Millan. New York. 1975).
- T. T. Puck. *Studies of the Life Cycle of Mammalian Cells*. Cold Spring Harbor Symposia. Vol. 29. pp. 167-176. (1964).
- J. J. Yunis. *Human Chromosome Methodology*. 2<sup>nd</sup> edition. p. 186. (Academic Press. New York. 1974).
- M. H. Wolf and K. E. Schnewies. *The Multiplication of Herpes Simplex Virus in the Different Phases of the Cell Cycle of Hep-2 Cells*. Zbl. Bakt. Hyg., I. Abt. Org. A. 224. pp.145-155.
- E. Robbins. *Cell Synchronization in Methods in Enzymology*. S. P. Colowick, N. O. Kaplan Ed. Vol XXXII Biomembranes Part B. p. 596. (1974).
- W. Nulisch. *Allgemeine Botanik Goerg Thieme Verlag*. Stuttgart. 6<sup>th</sup> edition. p. 328. (1971).
- H. E. Street. *Single Cell Clones in Street H. Plant Tissue and Cell Culture*. P. 203. (Blackwell Oxford. 1973).

FUJIFILM Irvine Scientific, Inc.  
2511 Daimler Street, Santa Ana, California 92705 USA  
Telephone: 1 949 261 7800 • 1 800 437 5706 Fax: 1 949 261 6522 • www.irvinesci.com

© 2019 FUJIFILM Irvine Scientific, Inc. All rights reserved. FUJIFILM Irvine Scientific, and the FUJIFILM Irvine Scientific logo, are trademarks of FUJIFILM Irvine Scientific, Inc. in various jurisdictions. All other trademarks are the property of their respective owners  
PN 40129 Rev.9

#### ENGLISH

##### INDICATION FOR USE

Colcemid Solution is intended for *in vitro* diagnostic use.

##### DEVICE DESCRIPTION

Colcemid Solution, is a ready-to-use solution with a concentration of 10 µg/mL of N-deacetyl-N-methyl colchicine (C<sub>21</sub>H<sub>25</sub>NO<sub>3</sub>) in Hank's Balanced Salt Solution. Colcemid is used to arrest mitosis in plant and animal cells by preventing spindle formation during metaphase of the cell cycle.<sup>1</sup> This effectively allows for the optimal yield of metaphases through enabling synchronization of cell cycles to facilitate cytogenetic analysis. Colcemid Solution is at least 30-40 times less toxic than colchicine.<sup>2</sup>

##### QUALITY ASSURANCE

A Safety Data Sheet and lot-specific Certificate of Analysis are available upon request.

##### SUGGESTED USES

- Chromosome Analysis – amniotic cells after initial cultivation.<sup>3</sup>  
Colcemid Solution arrests mitotic cultured cells in metaphase. Karyotyping for chromosome analysis is greatly facilitated when large numbers of metaphase cells are present in culture.
- Cell Synchronization  
Colcemid Solution aids in determining the mitotic activity index for cell synchronization studies.<sup>4</sup> Through investigations, Colcemid Solution has been shown to inhibit HeLa and CHO cells at concentrations of 0.05 – 0.5 µg/mL.<sup>5</sup>
- Polyploid Plants and Asexual Reproduction  
Colcemid Solution produces an increase in the chromosome count. New plant species can be produced in this manner under suitable conditions.<sup>6</sup> In asexual plant reproduction, the chromosome complement can be doubled under certain conditions.<sup>7</sup>

##### STORAGE INSTRUCTIONS AND STABILITY

Store tightly closed container at 2 – 8° C, protected from light.

When stored as directed, Colcemid is stable until the expiration date shown on the bottle label.

##### PRECAUTIONS AND WARNINGS

Colcemid Solution is mutagenic, tumorigenic, embryotoxic and teratogenic.

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