

Anti-Clumping Supplement

Catalog #	Product	Size
91150	Anti-Clumping Supplement	50 mL

Intended Use

This product is for research or further manufacturing use only.

Product Description

Anti-Clumping Supplement is designed to reduce cell aggregation for various cell types in suspension culture. This supplement is animal component and enzyme-free. This supplement is not intended to dissociate adherent cells.

Quality Assurance

All quality control test results are reported on a lot specific Certificate of Analysis which is available upon request.

Storage Instructions and Stability

Handle using aseptic techniques to avoid contamination. Store at room temperature or 2-8°C. Do not use any bottle of medium that shows evidence of particulate matter or cloudiness.

Precautions

The safety and efficacy of this product in diagnostic or other clinical uses has not been established. Please refer to the Safety Data Sheet for information regarding hazards and safe handling practices.

Directions for Use

Anti-Clumping Supplement is designed for use at a 1:200 – 1:1000 dilution (1 – 5mL/L) range. Typically used at 2mL/L. Supplement additional amount if necessary.

Note: Anti-Clumping Supplement will completely inhibit transfection via cationic liposomes or cationic polymers such as polyethylenimine (PEI). Cells must be spun down and re-suspended in media without Anti-Clumping Supplement before transfection.

2mL/L of Anti-Clumping Supplement added to HEK293 cultures

Without supplement



With supplement



Technical Support

CONTACT US

For more information or assistance contact Customer Service at:

- Email: fisitmrequest@fujifilm.com
- Direct line: +1 800 577 6097

WEBSITE RESOURCES

Visit the website at www.irvinesci.com for technical resources and information including:

- Safety Data Sheets (SDS)
- Certificate of Analysis (CoA) (when available)
- FAQs
- Product literature
- Complete list of offices and contact information by country

FUJIFILM Irvine Scientific, Inc.

2511 Daimler Street, Santa Ana, California 92705-5588 USA

Telephone: 1 949 261 7800 • 1 800 437 5706

Fax: 1 949 261 6522 • www.irvinesci.com

© 2020 FUJIFILM Irvine Scientific, Inc. All rights reserved. FUJIFILM Irvine Scientific and its logo are registered trademarks of FUJIFILM Irvine Scientific, Inc. in various jurisdictions. P/N 41062, Rev.01